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Downloads

[Overview](#)

Bead Lot Files for BD Cytometer Setup & Tracking Beads

Each lot of BD™ Cytometer Setup & Tracking beads has an associated bead lot file that needs to be downloaded from this website and then imported into BD FACSDiva™ software v6.0 or later.

The available bead lot files can be downloaded from the [BD FACSDiva™ Support](#) page.

Download Instructions

To properly download and install the appropriate bead lot file, please follow these instructions. (You may wish to print them.)

1. In the **Bead Lot Files** page linked above, click the link that matches the bead lot number on your vial of BD Cytometer Setup and Tracking beads. This will begin the download process.
2. If you are downloading directly to the cytometer workstation, save the zipped bead lot file to **C:\Program Files\BD FACSDiva Software\CST\Bead Lot**.

Otherwise, download the bead lot file to any location and manually transfer it via network, flash drive, or other method to: **C:\Program Files\BD FACSDiva Software\CST\Bead Lot**.
3. After placing the bead lot file in the proper folder, unzip it and start BD FACSDiva software on the cytometer workstation.
4. Within BD FACSDiva software, select **Cytometer > CST > Tools > Bead Lots**.
5. In the **Bead Lot** window, click the **Import** button.
6. Highlight the appropriate bead lot file and click **Open**.
7. Close the **Bead Lot** window.

Key Resources

[CS&T Bead Lot Files](#)

[Robust Statistics in BD FACSDiva™ 6.0 Software](#)

[BD™ CS&T Beads Data Sheet](#)

[All Resources >](#)

Related Products

[BD FACSCanto™ II Flow Cytometer](#)

More Information

[Ask BD](#)

[Sign Up for Email Updates](#)

Help

[Software Support](#)

[Technical Resources](#)

Support

[Overview](#) [Ordering](#) [Technical Support](#)

BD FACSDiva Software

Resources

Application Notes

[Bead Lot Files](#)

Product Information Sheets

[Quick Reference Guides](#)

Training

[Webinars](#)

Application Notes

• Qr and Br in BD FACSDiva™ v6 Software

• Establishing Optimum Baseline PMT Gains to Maximize Resolution on BD Biosciences Digital Flow Cytometers

• BD™ CompBeads plus BD FACSDiva™ Software: A Setup System

• BD FACSDiva™ Option White Paper

• Robust Statistics in BD FACSDiva™ 6.0 Software

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C:/Program Files/BD FACSDiva Software/CST/Beads Lot